

BarbLock™

Ultra-Secure Tubing Retainer

With 360° Radial Crimp & Seal



BarbLock on Coupling Insert by Colder Products



BarbLock on Coupling Body by Colder Products



BarbLock on 3/4" Mini Sanitary Fitting by Saint-Gobain Performance Plastics



BarbLock on 1.5" Sanitary Fitting by Saint-Gobain Performance Plastics



Cable Tie Replacement

NEW

Outperforms Cable Ties... Up To 7-1*


BarbLock Ultra-Secure Tubing Retainer Provides a High Performance Alternative to Cable Ties

- 360° Radial Crimp behind barb provides up to 7 times the pull-off resistance of cable ties
- 360° Compression Seal over the barb offers leak protection not found with cable ties
- Easy-to-assemble 2 piece design produces consistent retention and seal
- Materials used in the manufacture of BarbLock meet USP Class VI requirements
- Can be sterilized by radiation, steam or ethylene oxide
- Sizes available for most flexible tubing up to 3/4" outside diameter
- Recommended for use with Silicone, C-Flex®, Pharmed®, and PVC tubing products

Custom Sizes & Colors Available

- BarbLock Ultra-Secure Tubing Retainer can be sized to fit your existing barbed fitting
- Available for 1/8" to 1" hose barb
- Custom sleeve colors available

Contact your BarbLock representative for a Design Consultation.

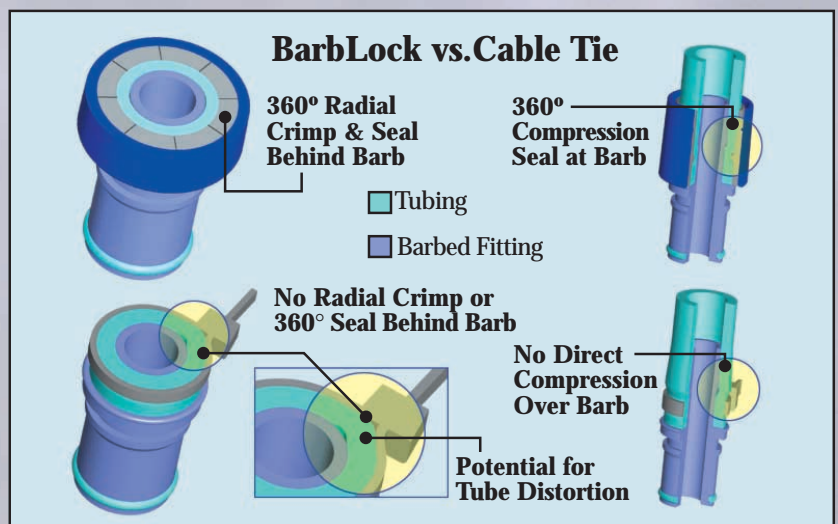


BarbLock Pump & Transfer Hose Assemblies




Clean room assembly and gamma sterilization available

BarbLock Hose Assemblies can be manufactured to your specifications for use as a peristaltic pump line or general fluid transfer line.


BarbLock Assembly

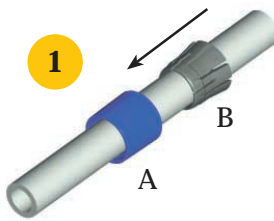


1
Install Sleeve (A) then Collet (B) over tube leading with ends marked "TOP".

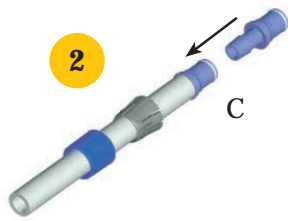

2
Install barbed end of fitting (C) into tube.

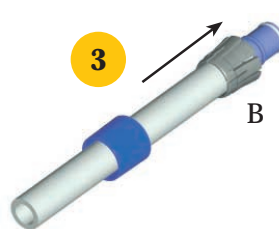


3
Lower Collet (B) over barb aligning base of Collet with end of tube.

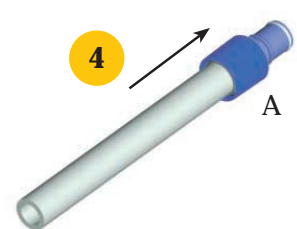


4
Using BarbLock Assembly Tool lower Sleeve (A) over Collet until locked.


BarbLock Performance vs. Cable Tie

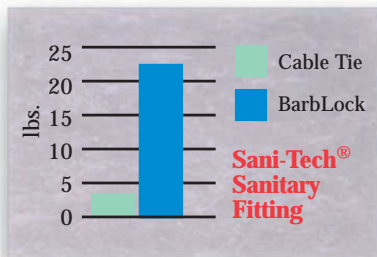

Pull-Off Strength with Silicone Tubing using Sani-Tech® 1.5" Sanitary Fitting

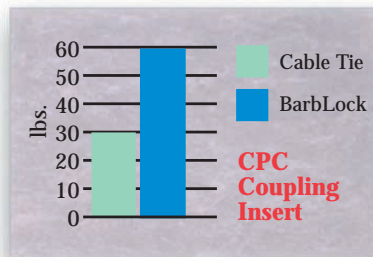

Pull-Off Strength with Silicone Tubing using Colder Products Coupling Insert

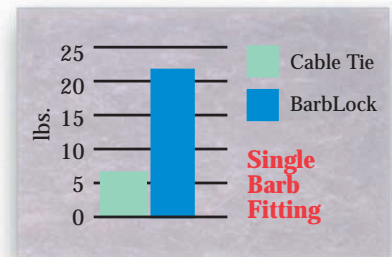

Pull-Off Strength with Silicone Tubing using Single Tier Polyethylene Barb

Distributed By:

Infiltec GmbH - Applied Critical Fluids
Alte Rheinhäuser Straße 8, 67346 Speyer
Germany

Tel. 0049/6232/24024
Fax 0049/6232/24025
www.infiltec.de - wz.infiltec@t-online.de



* 7-1 pull-off results attained using 1/2" ID x 3/4" OD 50 durometer silicone tubing and Sani-Tech 1.5" sanitary polypropylene fitting.

All products not manufactured by BarbLock Corporation carry the original manufacturer's warranty. Copies of original manufacturer warranties are available upon request. BarbLock Corporation makes no warranty, express or implied, with respect to products not manufactured by BarbLock Corporation. All products manufactured by BarbLock Corporation, are warranted to be free from defects at the time of shipment from BarbLock Corporation for a period of 6 months from the date of first sale. If a product is found to be defective by us, BarbLock Corporation only obligation will be to repair or replace the defective product, or at our option, refund the purchase price. User assumes all other risks, including risk of injury, loss or damage, whether direct, incidental or consequential, arising out of use, misuse, or inability to use any product supplied by BarbLock Corporation. BarbLock Corporation limited warranty is expressly in lieu of all other warranties, express or implied, including all implied warranties of merchantability and fitness for a particular purpose.

BarbLock Corporation, assumes no obligations or liability for any information or advice furnished by it, or for any results obtained with respect to these products. All such information and advice is given and accepted at buyers risk.

BarbLock is a trademark of TBM Corporation. C-Flex is a trademark of Consolidated Polymer Technologies. Sani-Tech is a trademark of Saint-Gobain Performance Plastics